

RECOMMENDED CLEANING PROCEDURE

Hettich Centrifuge – *Rotors and Accessories*

1. Clean with soap & water or a mild cleaning agent after every use (or at least once a week)
 2. If parts are contaminated by pathogenic or radioactive material we recommend:
 - For disinfection: *Helipur® H plus N* from *B. Braun Melsungen*
 - For the removal of radioactive material: *Decon neutracon®* from *Decon Laboratories Limited*
 3. Dry immediately after cleaning
 4. Lightly grease aluminium parts after drying using acid-free grease, e.g. vaseline
 5. With aerosol-sealed rotors and bio safety systems:
 - *Clean & check sealing rings regularly (weekly)*
 6. Rotor and Motor Shaft (clean once a month):
 - *Disassemble the rotor and clean*
 - *Lightly grease the motor shaft*
 7. Checked all parts on a monthly basis for corrosion damage
- ! Rotors and attachments may no longer be used upon indication of wear and tear or corrosion !***